

Instruction sheet


1. Product introduction

This product is of stepless speed to convey liquid by peristaltic pump, and widely used in medical, agriculture, chemical and so on.

2. Product list

Machine 1pcs, DC adaptor (24V) 1pcs, tube 1 meter

3. Product image

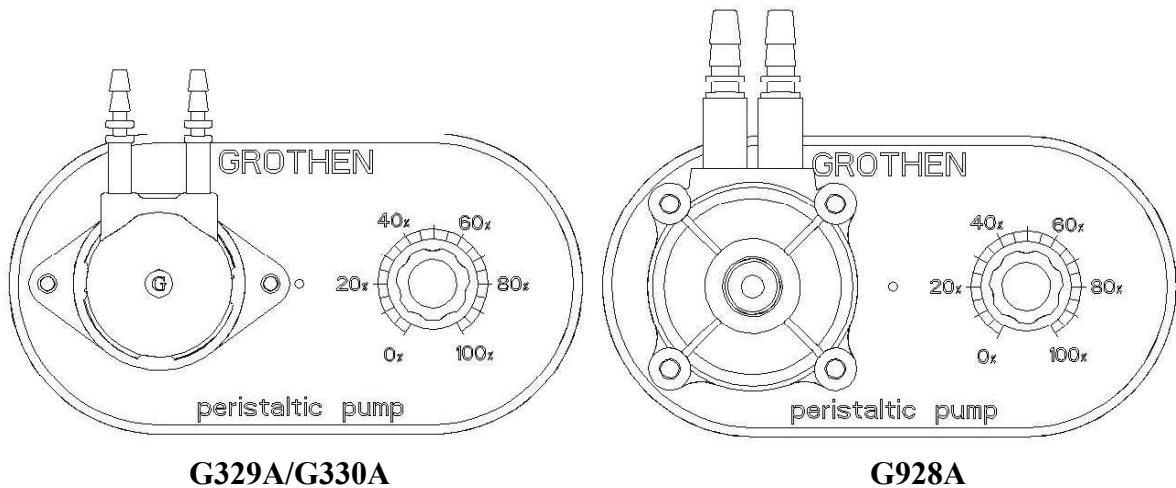

4. Product parameter

Model NO.	Pump head	Tube size : ID*OD mm	Tube Chemical Resistance	Tube life	Flowrate
G329A	1	3*5 food grade silicone tube	10% Acid / alkali	200H	15~95ml/min
G330A	1	2*4 pharmed bpt tube	30% Acid / alkali	1000H	6~60ml/min
G928A	1	6.4*9.6 food grade silicone tube	10% Acid / alkali	200H	114~500ml/min

5. Product use

1. Cut the tubes into two pieces and connect to the inlet and outlet of pump respectively.
2. Plug in the 24V adaptor.
3. Rotate the knob clockwise, and the pump works with speed bigger and with LED lighter gradually.
4. After use, rotate knob counter-clockwise to end, and then pump stops and LED goes out.